

Run-time error '75' or '339' FathZip.DDL in some of Vista OS

Cause:

- Download current version of DinoCapture (or V2.9.0.0 after)

If DinoCapture Version prior to V2.9.0.0
For Vista OS DinoCapture program required to set “Run as Administrator”
If you have any such problems, it is related to your Vista authorization issue.

Solution A:

- Right Click DinoCapture icon that is on the Desktop > select “Properties” > click on “Compatibility” tab > check “Run this program as an administrator” > then press “OK”

Solution B:

- When you run DinoCapture program for the first time, please don't double click on the DinoCapture icon. Instead, right click on the DinoCapture icon, then the optional manual will pop up and you should select "Run as administrator" Then confirm OK, the program will start to run. This is a one time process only, after that your computer will recognize that Run DinoCapture is authorized by the administrator. You will be able to double click DinoCapture for future use. If required to set “Run as Administrator” each time you start DinoCapture, please use solution A.

