

## Unable to save pictures / get an error message after taking a picture.

### Cause:

- ❏ Possible corrupted picture in “My documents”.

### Solution A:

- ❏ Delete the last picture taken with DinoCapture located in:  
(My documents > Digital Microscope > Picture)

### Solution B:

- ❏ Move or delete the pictures from "Digital Microscope" folder located in My documents (My documents > Digital Microscope > Picture) then start DinoCapture again and save images again to see if it works.

### Solution C:

- ❏ Another possible solution is to directly install DinoCapture on C:/ drive (eg. C:/Dino)

