

# ERROR CODE - RUNTIME ERROR 75

## Path/File access error

### Cause:



#### Vista OS

Program does not have rights or access to a file. For Vista OS required to set “Run as Administrator”

### Solution:



Right Click DinoCapture icon that is on the Desktop > select “Properties” > click on “Compatibility” tab > check “Run this program as an administrator” > then press “OK”



#### XP OS

Often this is caused when a program is trying to access a network file it doesn't have proper access to either because of network privileges or something is blocking the program. This issue can also be caused when the file is being used by another program or is read-only.

<http://www.computerhope.com/issues/ch000380.htm>

