

ERROR CODE - RUNTIME ERROR 50003

Unexpected error

Cause:

- Confusion or outdated Windows Visual basic or OCX controls

Solution:

- Try uninstalling DinoCapture and Dino-Lite drivers. Then close all open programs and reinstall our software again.

Source:

- http://www.rebrandsoftware.com/showsupport.asp?article_id=46
- Another solution is running windows update and installing both Critical and Optional software updates which might solve the issue.
- After updating your PC, try re-installing the DinoCapture software Refer-ences:
- <http://support.microsoft.com/kb/q194414/>
- <http://www.developersdex.com/vb/message.asp?p=640&r=5752013&Page=1>
- http://www.rebrandsoftware.com/showsupport.asp?article_id=46

