

Install DinoLite drivers each time after you reboot the computer in a network environment.

Solution:

- Have the Network Administrator install then run and close the program for the first time or at least one time.
- To Check if DinoCapture is already installed in your computer go to: Run
(type) REGEDIT
- Find the file located in-
Hkey_Local_machine
System
CurrentControlSet
Enum
USB
Vid_a168&Pid.....
- If the device is plugged in. It should show VID_A168*****
if unplugged, it should still be there.
- If the DinoLite is plugged in and it is still not working, then delete this file and reinstall it again.
It is possible that the file is corrupted.

