

The view window is larger than normal on DinoCapture (view window partially covers the “video” tab)

Solution A:

- Download current version of DinoCapture (or V2.9.0.0 after)

If DinoCapture Version under V2.9.0.0

Solution B:

- Please check your Display property >Setting >Advance Display
DPI 96 or 120? Please use 96 DPI.

Solution C:

- If reduced DPI to 96 but prefer larger font's you can enlarge by manipulating display properties:
Display property > “Appearance” tab > Font size: Extra Large Fonts
And to enlarge icons
Display property > “Appearance” tab > “Effects” button > check “Use Large Icons”

