

Windows 64 bit OS Compatibility issues

Solution A:

- Download current version of DinoCapture (or V2.9.0.0 after)

Solution B:

- You can still use 64 bit software that you currently use in conjunction with the Dino-Lite drivers, such as webcam software or video recording software. If you do not have any related software, you can try third party free or open source software that is available online such as and not limited to:

Dorgem <http://dorgem.sourceforge.net/>

or

AMCAP <http://amcap.en.softonic.com/>

- A note that since these are third party software, some features advertised might not be functional other than the basic use of the Dino-Lite Microscopes such as viewing and taking pictures.

