

How to remove DinoCapture program from XP or Vista

How to remove DinoCapture program from Windows VISTA

Unplug the Dino-Lite USB Cable then go to “Start” located on the desktop > control panel > uninstall a program > Look for Dino Lite Plus and Pro > then select “uninstall”. Then go back to “start” menu on desktop > go to “All Programs” > then look for “Dino Digital Microscope” folder and expand the folder > select “Uninstall DinoCapture” If you have taken video or picture before go to the “start” menu and then the “Documents” folder. Back up the pictures or delete the “Digital Microscope” folder.

How to remove DinoCapture program from Windows XP

Unplug the Dino-Lite USB Cable then uninstall the Dino-Lite drivers through the control panel:
Start~> Control panel~>add or remove programs. Then delete DinoCapture by going through the “Start” menu~> All Programs~>Choose “Dino Digital Microscope” program~> Scroll to “Uninstall Driver” Now back up the pictures and video that you want to keep into another folder and delete the folder called “Digital Microscope” located in “My Documents” folder.

